INSTRUCTIONS FOR USE

ACTIVADISC® Depth Filter Modules

Installation of ACTIVADISC® depth filter modules

Before use, check the filter modules conditions including gaskets. Defective filter modules should not be used. The modules must be inserted centrally according to their adapter and shall not get in contact with the housing at any point.

Rinsing the filter modules

- Use 50 L/m² of a process-compatible solution (e. g. WFI, ethanol)
- The rinsing flow rate shall be approximately 350 L/m²×h
- Rinse in filtration direction (forward flow)

If the rinsing is not desired or tolerated (insoluble products such as high percentage distillates or oils), the product to be filtered might be first recirculated to remove possible small fibers and/or diatomite particles.

Sterilization

- Fill the system with water, slightly open all outlet valves and increase the temperature gradually
- Heat up the filter until the required sterilization temperature of 85 °C is reached at all outlet valves
- Maintain temperature for minimum 20 minutes
- Turn off the hot water supply
- Cool down gradually the filter to room temperature with cold sterile water while recirculating it into the system

To save energy, it is recommended in e) and f) to heat up during recirculation.

NOTE

Sterilization is not fully completed until all parts that come into contact with the product have been sterilized. For alternative sterilisation methods, please contact your FILTROX representative.

Filtration

- Slowly start filling the filter by opening the inlet valve and all venting valves (ensure complete venting), the outlet valve should be closed
- Once properly vented, open the outlet valve, close the venting valves and start the filtration (if required, the initial filtered volume may be separated from filtrate to avoid any dilution with rinsing water)
- Adjust the flow rate to 150 L/m²×h or according to the process needs
- Ensure continuous and uniform flow rate
- Avoid any abrupt changes in filtration flow, pressure shocks, gas bubbles as well as stop and go

NOTE

Observe the maximum recommended flow rate, operating pressure and temperatures. Further information can be found in the corresponding technical data sheet.

Regeneration

Once the filters are saturated and the filtration-treatment is no longer in accordance with requirements, the ACTIVADISC® depth filter modules are to be considered exhausted and cannot be regenerated.

Replacement of filter modules

The ACTIVADISC® depth filter modules must be replaced as soon as the filtration-treatment requirements are not fulfilled. The filtered volumes vary depending on products, process target, flow rate as well as ACTIVASORB® grade.

Note on disposal

The respective official regulations for disposal must be followed, depending on the filtered product. Uncontaminated modules can be disposed of as non-hazardous waste. Please refer to your local environmental authority for more information.

Shelf life and storage

The filter modules should be stored in their original packaging in a dry, odor-free and well-ventilated area. Temperature fluctuations can lead to condensation formation and should therefore be avoided. The filter modules are intended for immediate use. Filter modules for food & beverages should be used within 60 months from date of manufacture. Filter modules for life science applications should be used within 36 months from date of manufacture. Further information is given in the corresponding technical data sheet for filter media.

Technical Support

Contact your local FILTROX representative or our experts at applications@filtrox.com. We will be happy to help you to find the best possible filtration solution for your product.

Remarks

All information is based on current knowledge and does not claim to be complete. No liabilities can be derived from this information. FILTROX reserves the right to make changes in the course of technical improvement.

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